


Gravimetric Determination of NaHCO₃ in a Mixture

Final Exercise - 100 points

100 / 800

1

4. What is the stoichiometry of the reaction?

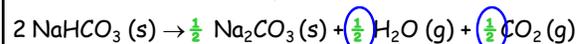

$$n_{\text{CO}_2} = n_{\text{H}_2\text{O}} = \frac{1}{2} n_{\text{NaHCO}_3}$$

$$\text{but, weight loss} = n_{\text{CO}_2} * 44.01 + n_{\text{H}_2\text{O}} * 18.02$$

$$\text{so, weight loss} = \frac{1}{2} n_{\text{NaHCO}_3} * (44.01 + 18.02)$$

$$n_{\text{NaHCO}_3} = 2 * (\text{weight loss}) / 62.03 = (\text{weight loss}) / 31.02$$

$$5. w_{\text{NaHCO}_3} = n_{\text{NaHCO}_3} * 84.01 \text{ g/mol}$$

Molar Mass of NaHCO₃

6. Compute Percent Composition of Sample

$$\text{Pct}_{\text{NaHCO}_3} = 100 * w_{\text{NaHCO}_3} / w_{\text{Sample}}$$

4

Purpose:

Determine the percent of NaHCO₃ in a mixture



Concepts:

Thermal Decomposition
Constant Weight

Complete Reaction
Limiting Reagents

Techniques:

Weighing by Difference
Thermal Decomposition



Apparatus:

Analytical Balance
Crucible

Hot Plate
Crucible Tongs



Gravimetry: An analytical technique based primarily on weighing

2

Summary: Pct NaHCO₃ is proportional to the weight loss

Showed that:

$$n_{\text{NaHCO}_3} = (\text{weight loss}) / 31.01$$

$$w_{\text{NaHCO}_3} = n_{\text{NaHCO}_3} * \text{MM} = ((\text{weight loss}) / 31.01) * 84.01$$

$$= 2.709 * (\text{weight loss})$$

$$\text{Pct}_{\text{NaHCO}_3} = 100 * w_{\text{NaHCO}_3} / w_{\text{Sample}}$$

$$= 270.9 * (\text{weight loss}) / w_{\text{Sample}}$$

$$w_{\text{NaHCO}_3} = (\text{weight loss}) / 0.3691$$

What weight would be lost by 1.0000 g of pure NaHCO₃?

$$\text{weight loss} = \text{Pct}_{\text{NaHCO}_3} * w_{\text{Sample}} / 270.9$$


$$\text{weight loss} = 0.3691 \text{ g per g of NaHCO}_3$$

5

What we do.

1. Weigh Sample, w_{Sample} .

2. Decompose the sample by heating

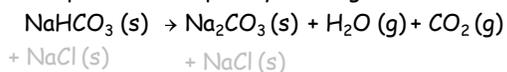

3. Weigh product

$$\text{weight loss} = w_{\text{CO}_2} + w_{\text{H}_2\text{O}}$$

$$\text{weight loss} = n_{\text{CO}_2} * 44.01 + n_{\text{H}_2\text{O}} * 18.02$$

weight = mol X Mol Mass

Molar Mass of CO₂

Molar Mass of H₂O³

What Can Go Wrong?

The accuracy and precision of the result depend on:

- care with which weighings are done
 - weigh unknowns by difference
 - weigh only cool samples
 - don't touch the crucible
- preventing loss of sample by spilling
 - don't spill
- care with which decomposition is conducted
 - make sure reaction is complete
- calculation or transcription errors
 - don't complete data sheet before end of lab
 - check your arithmetic and data entries

6

This exercise depends heavily on the proper use of the analytical balance!

- try to use same balance throughout a run

DO NOT TINKER WITH BALANCE SETTINGS!

?? Standard Penny ??

Weigh a coin before anything else

Check the weight of the "standard" coin before each subsequent weighing.

7

How Do We Know When the Reaction is Complete?

When all the H₂O and CO₂ have been lost, the sample will cease to lose weight. So we weigh and re-heat.

If the reaction is complete, two successive weighings before and after heating should be the same

NOTE
±

How should we define "the same" (δ)?

NOT
0.5 mg!

Our operational definition is that they are the same if they differ by less than ± 5 mg. I.e., $|w_2 - w_1| < 0.0050$ g

e.g.	16.5386	w ₂	16.5294	w ₃	16.5242	w ₃	16.5242
w ₂	<u>16.5294</u>	w ₃	<u>16.5242</u>	w ₄	<u>16.5214</u>	w ₄	<u>16.5276</u>
	- 0.0092		- 0.0052		- 0.0028		+ 0.0034
	- 9 X mg		- 5 X mg		- 2.8 ✓ mg		+ 3.4 ✓ mg

Weighing by Difference

Most errors in weighing are due to loss of material in the transfer from one container to another!

How do we minimize this problem?

- Minimize the number of transfers
- Don't use intermediate containers or devices



8

How do the Various Substances behave at High Temperature?

NaHCO₃ decomposes rapidly at T > 200 °C forming Na₂CO₃

NaCl melts at 800 °C, but does not decompose.

Na₂CO₃ (the product) melts at 850 °C AND begins to decompose to CO₂ + Na₂O at that temperature.

Would result in additional weight loss

850 °C is bright red heat



But make sure it is at least 200 °C

The peak temperature on a hot plate is well below 800 °C.

11

Weighing by Difference

- Process:
 - weigh sample container,
 - transfer sample *directly* into final container by tapping
 - reweigh original sample container

- Repeat until
 - Difference between initial and final weights of container is the desired sample weight

You are NOT "weighing by difference" if you:

- use a spatula
- place heavy flask/beaker on balance pan
- use a watch glass or piece of paper
- record (only) weight of sample - not $w_f - w_i$

Web includes a slide show describing weighing by difference.

9

Reproducibility

Uncertainty in Weighing:

w_{sample} analytical balance (0.0002 / 1.0000) < 0.02%

w_{residue} analytical balance (0.0002 / 0.7000) < 0.02%

Our criterion for complete reaction is weights that differ by less than 5 mg. So, our weight loss should differ from the "true" weight loss by less than 5 mg.

Unknowns are 50 - 100% NaHCO₃. We weigh 1.5 - 2.0 g

A 2 g sample with 50% NaHCO₃ will contain 1 g of NaHCO₃ and the (*minimum*) weight loss will be 0.3691 g

5 mg constitutes less than a $100 \times 0.005 / 0.3691 = 1.4\%$ error in weight loss

Any larger error will be due to loss of sample or incomplete thermal decomposition.

12

Do the analysis twice,
and

Do Not Burn Yourself

Handle the crucible **ONLY** with crucible tongs!

13

DATA SHEET

Wt of crucible + sample	16.0755 g	
Wt of crucible	13.9842 g	399.5 mg
Wt of sample	2.0913 g	
Wt of crucible + residue –aft heat1	15.6760 g	19.7 mg
Wt of crucible + residue –aft heat2	15.6563 g	9.3 mg
Wt of crucible + residue –aft heat3	15.6470 g	3.4 mg
Wt of crucible + residue –aft heat4	15.6436 g	
Wt loss [16.0755 – 15.6436]	0.4319 g	
Wt NaHCO ₃ [0.4319 / 0.3691]	1.170 g	Wt loss per gram of NaHCO ₃
% NaHCO ₃ [100 X 1.170 / 2.0913]	55.95 %	

14

Review - Avg, Avg Deviation, Pct Deviation

Our first result is 55.95%. Suppose our second result is 56.42%

Average = $\frac{55.95 + 56.42}{2} = 56.19 \%$

Avg Dev = $\frac{|55.95 - 56.19| + |56.42 - 56.19|}{2}$

= $\frac{0.24 + 0.23}{2} = 0.24 \%$

Pct Dev = $\frac{100 \times 0.24}{56.19} = 0.43 \%$

15